


InertSustain™ Phenylhexyl

Physical Properties


● Silica	: New Evolved Surface (ES) Silica Gel	● Bonded Phase	: Phenylhexyl Groups
● Particle Size	: 3 μm, 5 μm	● End-capping	: Yes
● Surface Area	: 350 m ² /g	● Carbon Loading	: 9.0 %
● Pore Size	: 100 Å (10 nm)	● USP Code	: L11
● Pore Volume	: 0.85 mL/g	● pH Range	: 1.0 ~ 10.0

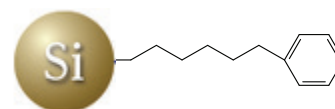


InertSustain Phenylhexyl columns are bonded with Phenylhexyl groups, which employs a Phenyl ring with a Hexyl (6- Carbon) linker and is Densely bonded to our newly developed Evolved Surface silica gel delivering complementary selectivity to straight Alkyl-chain columns, but with industry leading inertness, lot-to-lot reproducibility and low back pressure.

ODS column changed the separation pattern

Phenyl column performed a strong retention

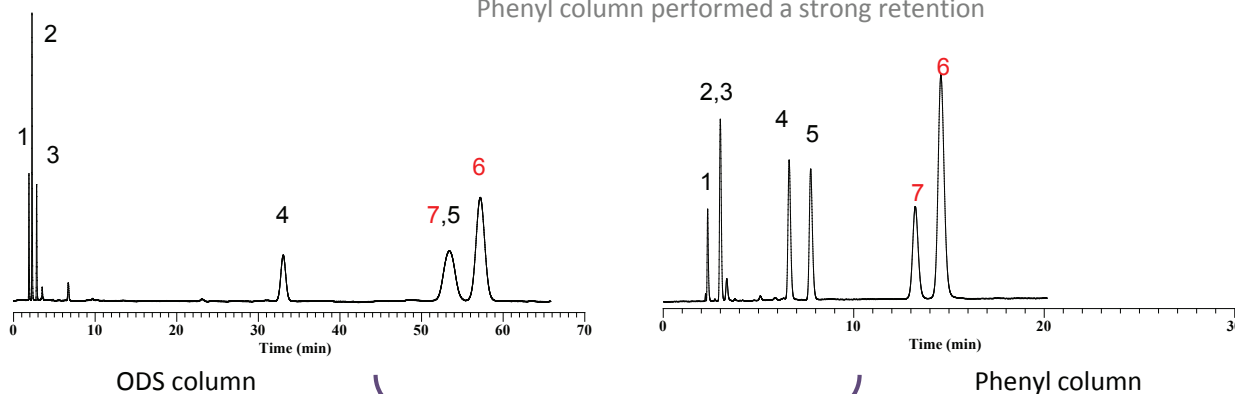


InertSustain Phenylhexyl

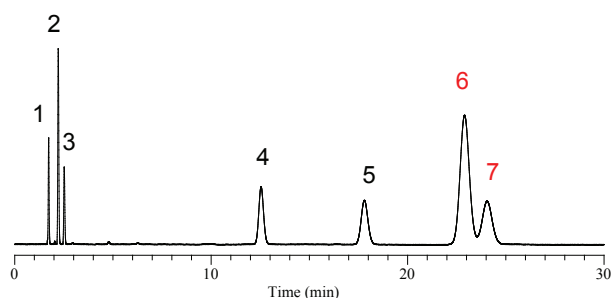


Conditions

Eluent	: A) CH ₃ OH, B) H ₂ O, A / B = 70/30, v / v
Flow Rate	: 1.0 mL/min
Col. Temp.	: 40 °C
Detection	: UV 254 nm
Sample	: 1. Uracil 2. Caffeine 3. Phenol 4. n-Butylbenzene 5. n-Amylbenzene 6. Triphenylene 7. o-Terphenyl

Analysis of Strong Acidic Compound

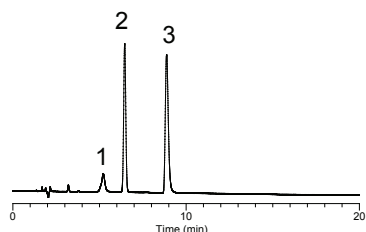


InertSustain Phenylhexyl

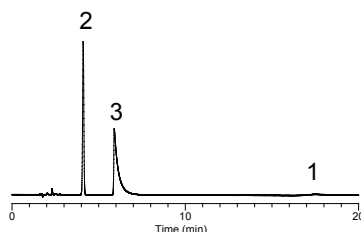

XBridge Phenyl

Conditions
 Column Size : 5 µm, 150 x 4.6 mm I.D.
 Eluent : A) CH₃CN
 B) 0.1% H₃PO₄
 A/B = 25/75, v/v
 Flow Rate : 1.0 mL/min
 Col. Temp. : 40 °C
 Detection : UV 254 nm
 Sample : 1. Brilliant Blue FCF
 2. Phenol
 3. Salicylic acid

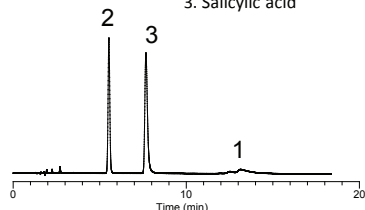

Luna Phenyl-Hexyl

Ordering Information:

HP Series Particle Size: 3 µm 50 Mpa(500 Bar)	Length \ I.D. (mm)	2.1	3.0	4.6
		30	5020-89209	5020-89215
	50	5020-89210	5020-89216	5020-89222
	75	5020-89211	5020-89217	5020-89223
	100	5020-89212	5020-89218	5020-89224
	150	5020-89213	5020-89219	5020-89225
	250	5020-89214	5020-89220	5020-89226

Particle Size : 3 µm	Length \ I.D. (mm)	1.0	1.5			
		30	5020-89160	5020-89166		
		50	5020-89161	5020-89167		
		75	5020-89162	5020-89168		
		100	5020-89163	5020-89169		
		150	5020-89164	5020-89170		
		250	5020-89165	5020-89171		
	Length \ I.D. (mm)	2.1	3.0	4.0	4.6	
		30	5020-89124	5020-89131	5020-89138	5020-89145
		50	5020-89125	5020-89132	5020-89139	5020-89146
		75	5020-89126	5020-89133	5020-89140	5020-89147
	100	5020-89127	5020-89134	5020-89141	5020-89148	
	150	5020-89128	5020-89135	5020-89142	5020-89149	
	250	5020-89129	5020-89136	5020-89143	5020-89150	

Particle Size : 5 µm	Length \ I.D. (mm)	1.0	1.5			
		30	5020-89038	5020-89044		
		50	5020-89039	5020-89045		
		75	5020-89040	5020-89046		
		100	5020-89041	5020-89047		
		150	5020-89042	5020-89048		
		250	5020-89043	5020-89049		
	Length \ I.D. (mm)	2.1	3.0	4.0	4.6	
		30	5020-89001	5020-89008	5020-89015	5020-89022
		50	5020-89002	5020-89009	5020-89016	5020-89023
		75	5020-89003	5020-89010	5020-89017	5020-89024
	100	5020-89004	5020-89011	5020-89018	5020-89025	
	150	5020-89005	5020-89012	5020-89019	5020-89026	
	250	5020-89006	5020-89013	5020-89020	5020-89027	

* End-fittings are 1/16" Waters-compatible

Worldwide Ordering Information

GL Sciences Inc. Japan
 22-1 Nishishinjuku 6-Chome
 Shinjuku-ku, Tokyo, 163-1130, Japan
 Phone: +81-3-5323-6620
 Fax: +81-3-5323-6621
 Web: www.glsciences.com
 Email: world@gl.co.jp

GL Sciences B.V.
 De Sleutel 9, 5652 AS Eindhoven
 The Netherlands
 Phone: +31 (0)40 254 95 31
 Orders: orders@glsciences.eu
 Web: www.glsciences.eu
 Email: info@glsciences.eu

GL Sciences Inc. USA
 4733 Torrance Blvd. Suite 255
 Torrance, CA 90503
 Phone: 310-265-4424
 Fax: 310-265-4425
 Web: www.glsciencesinc.com
 Email: info@glsciencesinc.com

Distributor:

