

FastRemover™ for Phospholipid

Rapid and Efficient Removal of Proteins and Phospholipids

The FastRemover for Phospholipid 96-well plate deliver a rapid and effective removal of proteins and phospholipids in plasma and serum samples without sacrificing the recovery of your target analytes.





Benefits

- Simple & easy protocol to remove proteins and phospholipids
- High sensitivity analysis is unaffected by contamination from plasticizers or other impurities found in other 96-well plates
- Removal of microparticle contaminants enables injection to LC/MS/MS directly from the collection plate
- Removes more than 90% of phospholipids resulting in eliminating ion-suppression
- Prolong HPLC/UHPLC column lifetime by removing proteins and phospholipids that can damage your column

Phospholipid Removal

The presence of phospholipids in plasma or serum samples is one of the major problems in LC/MS-(MS) analysis. Phospholipids can build up on your MS system and bleed off the HPLC/UHPLC column, causing ion suppression, shifts in retention time and peak shape and necessitating time consuming column and system maintenance. Use of FastRemover for Phospholipid 96-well plate will eliminate these effects and extend the lifetime of your HPLC/UHPLC column and deliver more predictable/accurate mass spectrometry results.


Easy Procedure by FastRemover for Phospholipid

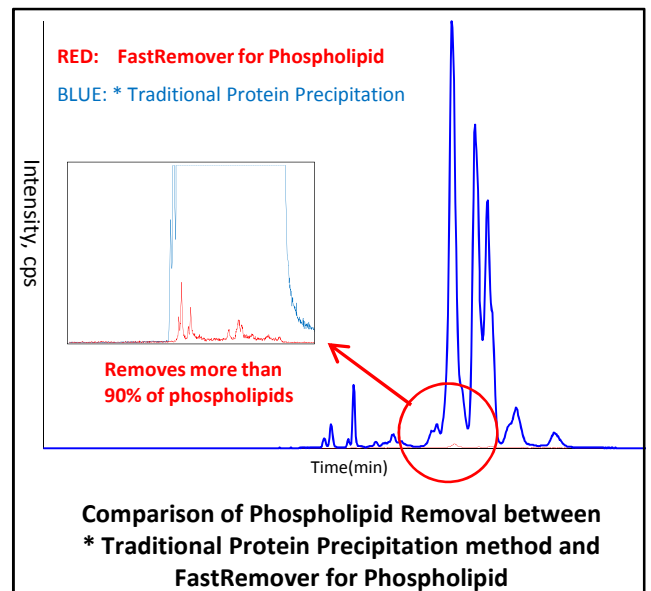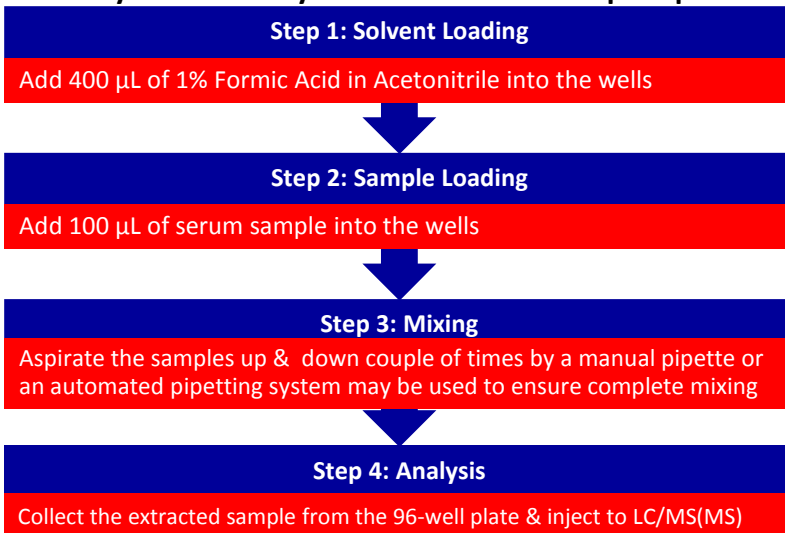

* Traditional Protein Precipitation method:
Adding acetonitrile to sample and collecting the supernatant layer


Extend HPLC/UHPLC Column Lifetime

Over the course of multiple injections, phospholipids build up and can lead to reduced column lifetime, showing increase in column back pressure, decrease in column sensitivity and efficiency. The figure on the right illustrates the removal efficiency of phospholipids, proteins and microparticles by FastRemover for Phospholipid.

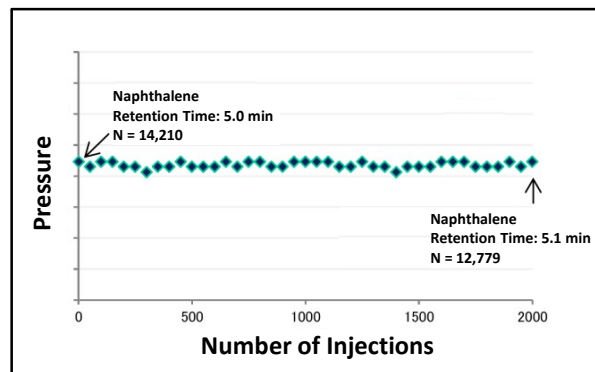

Industry Leading High Recovery for Bioanalysis

As shown below, the FastRemover for Phospholipid 96-well plate deliver a rapid and effective removal of proteins and phospholipids in plasma and serum samples without sacrificing the recovery of your target analytes.

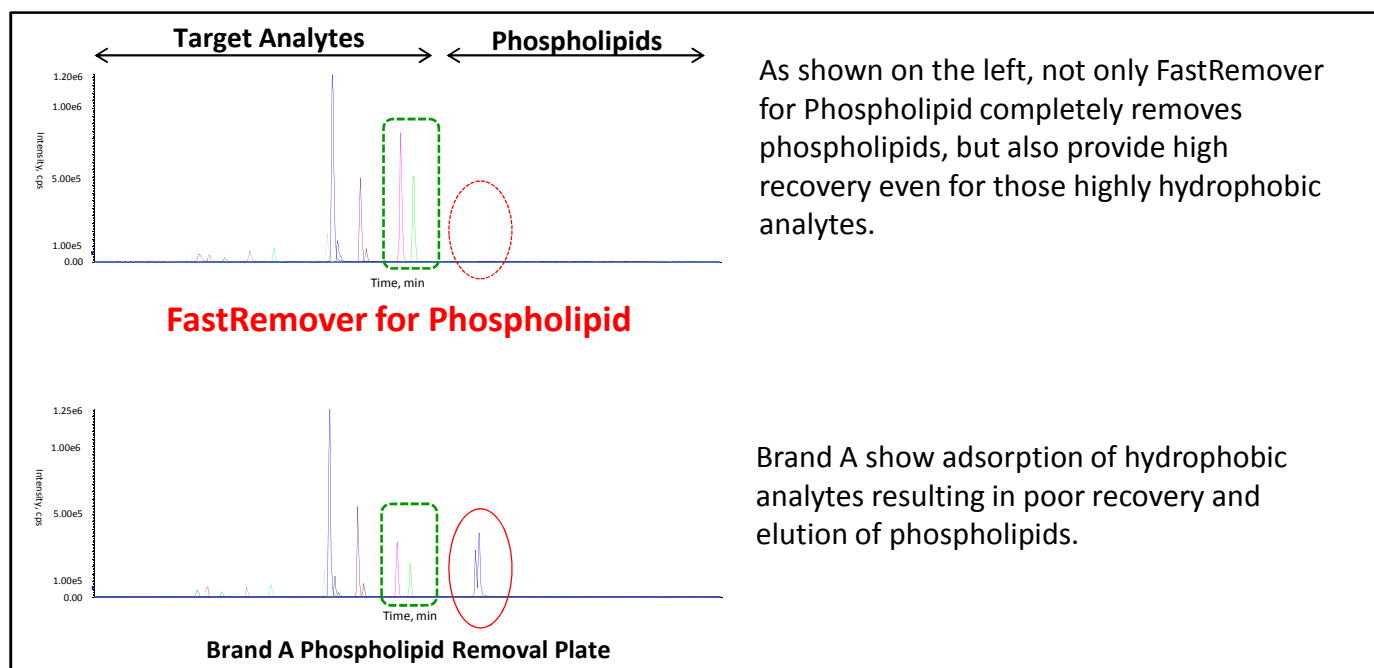

Ordering Information

Description	Qty/Pkg	Cat.No.
FastRemover for Phospholipid (0.2 μm)	1	7510-11021

Contact Information

GL Sciences, Inc. USA
 4733 Torrance Blvd. Suite 255
 Torrance, CA 90503
 Phone: 310-265-4424
 Fax: 310-265-4425
 www.glsciencesinc.com
 email: info@glsciencesinc.com

GL Sciences B.V.
 De Sleutel 9
 5652 AS Eindhoven
 The Netherlands
 Phone: +31 (0) 40 254 95 31
 www.glsciences.eu
 email: info@glsciences.eu

GL Sciences, Inc. Japan
 22-1 Nishishinjuku 6-Chome
 Shinjuku-ku, Tokyo, 163-1130, Japan
 Phone: +81-3-5323-6620
 Fax: +81-3-5323-6621
 www.glsciences.com
 email: world@glsc.co.jp

International Distributors
 Visit our Website at
www.glsciences.com/distributors